**DOUCHE DE SÉCURITÉ ANTIGEL
AXION® MSR EMERGENCY FREEZE PROTECTED SHOWER**

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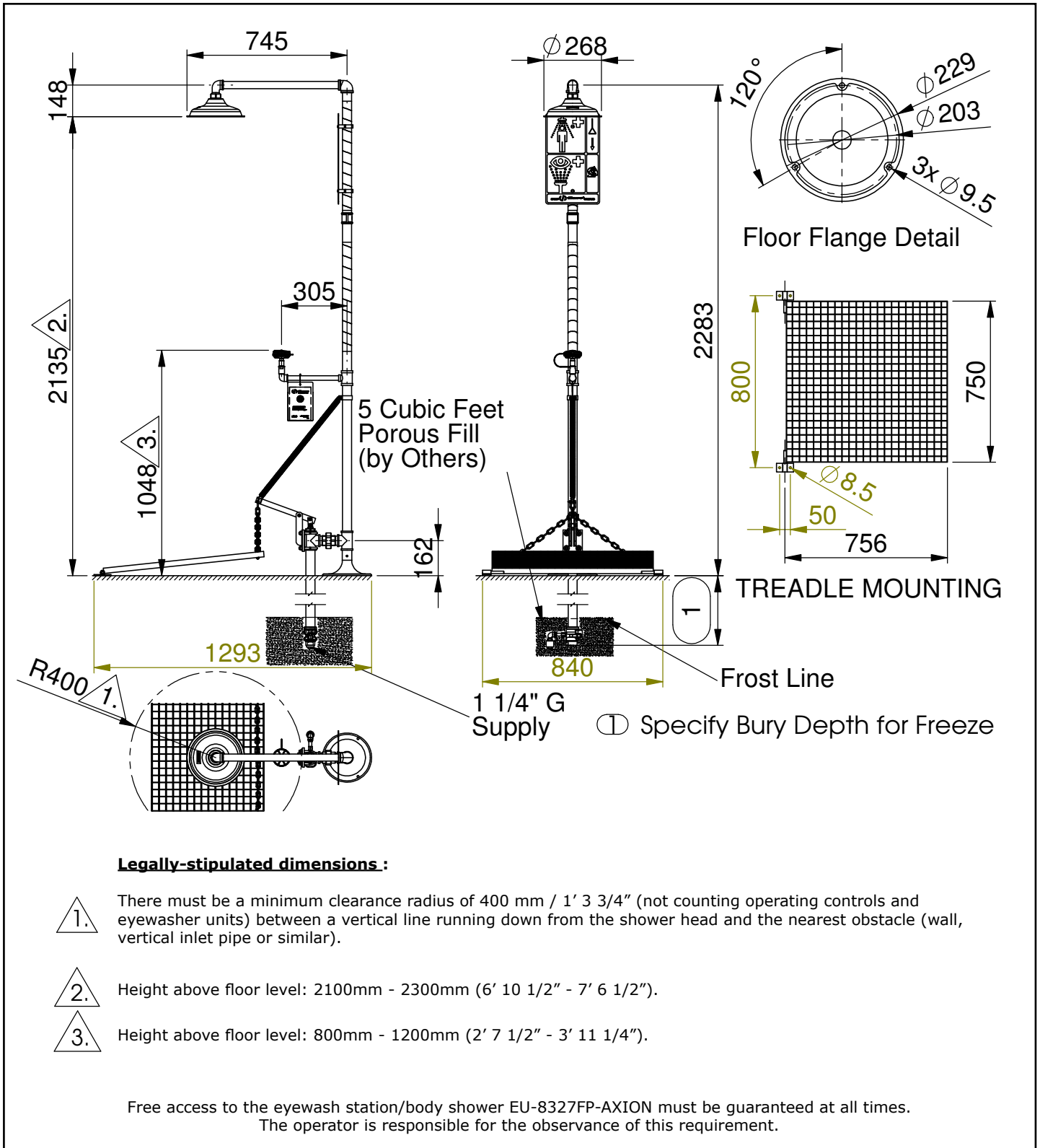
Model **EU-8327FP-AXION**, freeze protected combination shower and eye/face wash with AXION® MSR eye/face wash and showerhead, with a stainless steel grate.

Les dimensions peuvent varier +/- 12,7 mm (1/2")
These dimensions may vary +/- 1/2"

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Shower location :

The eyewash station/body shower EU-8327FP-AXION must be installed in the vicinity of potential danger zones. It must be highly visible and easily accessible.

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You are recommended to use a 1 1/4" G conduit, capable of withstanding pressures of 2 - 6BAR (30 - 90PSI). The use of an inlet filter designed to prevent the entry of suspended particles and other impurities is also recommended.

Water quality :

Eyewash stations/body shower require a supply of drinking water conforming to European standards or their equivalents in other countries.

Water temperature :

The required water temperature should be determined the user. The accepted ideal range is 15 - 35°C / 60 - 95°F (but please ensure conformity with local regulations).

Volume flow :

A minimum flow rate of 60 litres / 15,8 US gallons per minute must be maintained for the body shower. A minimum flow rate of 6 litres / 1,6 US gallons per minute must be maintained for the eyewash station. The eyewash station/body shower must be capable of delivering this amount of water for at least 15 minutes.

Connection to plumbing :

The inlet consists of a 1 1/4" G conduit with female-thread adapter.

Regular inspection :

The eyewash station/body shower EU-8327FP-AXION should be inspected at weekly intervals. The inspection tag should be dated and signed after each check.

TROUBLESHOOTING	
PROBLEM	REPAIR CHECKLIST
1. No flow	Check the main shutoff valve.
2. Insufficient flow of water	1. Check the line pressure. This should be at least 2 BAR / 30 PSI (dynamic). 2. Blocked flow, possibly due to the conduit not being correctly flushed out. Dismantle the flow-control element, and clean the filter insert.
3. Flow rate is not as specified	The flow conduit may be blocked (for remedy, see above). The eyewash station may not have been set up correctly.
4. Water does not discharge correctly	Check the building's main drainage outlet. Note that it must have adequate capacity to handle the entire drainage system.
5. Loose connections or fittings	Tighten any loose connection elements. The application of hot water sometimes makes it easier to put slightly-misaligned connection elements back into place. If the join is still leaking after 30 minutes, the assembly must be dismantled and resealed.

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